
🔸 Fix Me!

This page is not fully translated, yet. Please help completing the

translation.

(remove this paragraph once the translation is finished)

130mm Apochromat from Telescope-Service

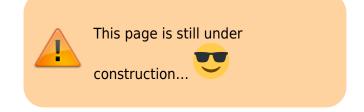

This page is about our TS-Optics-PHOTOLINE Apochromat. It has an aperture of 130mm and a focal length of 910mm. Consequently, the focal ratio is f/7 but can be reduced to f/5.2 with a reducer. It is equipped with a triplet air-spaced objective lens featuring an SFPL-53-ED glass element from Japan and a 3.7" focuser.

Eyepiece-side connections:

- 1.25″
- 2″
- M68x1 internal thread
- M92x1 internal thread

Available Accessories:

- TS-Optics 1.0x Flattener Gen 2 (TSFlat3)
- TS-Optics 0.74x Reducer (TSAPORED075)
- Baader Solar Filter for Optical Observation
- Baader Herschel Prism Mark II

Suitable Mounts:

Articles related to specific mounts are linked, and their setup is often described in combination with one of our other telescopes. However, the basic steps for assembling each mount are universal.

- Celestron CGE-Pro
- Celestron CGX-L
- Sky-Watcher EQ8-R Pro

Working Distance:

- 123mm from the 2" eyepiece holder
- can be extended to 223mm



Configuration

0.74x Reducer

Reducer + QHY600M with Filter Wheel and Off-Axis Guider

This setup is the standard for deep-sky observations as it provides the widest field of view. The reducer used is the TS-Optics 0.74x Reducer with:

- camera side M48x0.75 external thread with 55mm working distance
- camera side M76x1 external thread with 105.7mm working distance
- telescope side M88x1 external thread and M82x1 internal thread

The reducer reduces the focal ratio to f/5.2. The "drawback" of this setup is that the extension tube between the tube and the focuser must be removed, otherwise, achieving focus becomes impossible. This process requires care as the threads are delicate. For a comparison, see the following two images:



Correct setup without extension tube

Incorrect setup with extension tube

M68a → M62a Adapter

This setup requires the following adapters to be mounted in the order seen from the telescope:

TS-Optier 0.74x Reducer (TSAPORESOCI) M88 10mm etimeter (ute	Position	Adapter	Optical Path Length [mm]
	1	TS Optics M76 → M68 Adapter	20.7
ADD with Backward OUNGOD with Filter Miles and	2	M68 Fine-Tuning Ring (golden)	1
	3	M68a → M68i Conversion Ring	12
APO with Reducer, QHY600 with Filter Wheel and Off-Axis Guider, and all necessary adapters	4	M68 Fine-Tuning Ring (red)	0.5
The camera along with the filter wheel, filter,	5	M68 10mm Extension	10
and off-axis guider has a back focus of 56mm. The working distance behind the M76 thread of	6	M68 Fine-Tuning Ring (red)	0.5

7

The working distance behind the M76 thread of the reducer is 105.7mm. Thus, a distance of


5

49.7mm needs to be bridged. Simultaneously, a translation from M76 to M62 thread is required. The former is achieved with an M76 to M68 adapter from TS Optics, a conversion adapter from M68a to M68i, and an adapter from M68a to M62a. These three adapters have an optical path length of 37.7mm. The remaining 12mm are bridged by a 10mm M68 extension and three M68 fine-tuning rings with a total optical path length of 2mm.



Adapter combination as in the table (ascending from left to right)

It is important to note that before mounting the ¹ above adapter combination, the long M48 adapter, usually present on the reducer, must be unscrewed to expose the M76 thread.

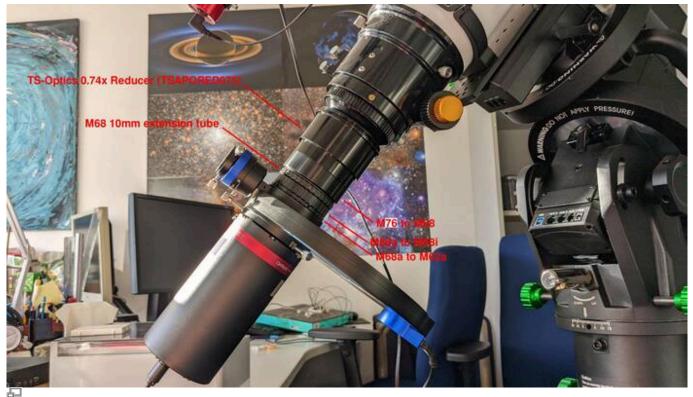

Reducer + QHY600M with UFC Filter Holder

Placeholder image (must be replaced)

The same principles apply here as mentioned above, with the adapter combination being slightly different:

Position	Adapter	Optical Path Length [mm]
	TS Optics M76 → M68 Adapter	20.7
2	M68 7.5mm Extension	7.5

Last update: 2024/02/19 en:ost:telescope:ts-apo https://polaris.astro.physik.uni-potsdam.de/wiki/doku.php?id=en:ost:telescope:ts-apo&rev=1708382738 22:45

ition	Adapter	Optical Path Length [mm]	
	M68a → M68i Conversion Ring	12	
4	UFC S70 / M68x1	1	
5	UFC Varilock	15-20	
6	UFC Tilter	9.75-10.5	
7	UFC Base	13	
8	Camera Adapter M54 → Camera Clamp	6	Placeholder image (to be replaced)

This adapter combination has an optical path length of 87.2 mm when the Varilock is set to 16.5-17.25 mm. This, combined with the QHY600M, which has a back focus of 17.5 mm, and the installed filters, results in a total working distance of 105.7 mm behind the M76 thread of the reducer.

QHY600M

The QHY600M can also be mounted on the APO without additional filter holders, filter wheels, etc. Only the QHY camera adapter with M54 internal thread and a 2" eyepiece holder from QHY with M54 external thread on one side and M84 internal thread on the other side are necessary. The M48 thread then fits onto the M84 thread of the reducer.



Position	Adoptor	Length [mm]	Adapter combination as in the table (ascending from left to right). Additionally, the adapter of
1	QHY Camera Adapter	5	the reducer with the M48 thread is mounted on
2	2" Eyepiece Holder	33	the right.

Together with the 17.5 mm back focus of the camera, this achieves the required working distance from the M48 of the reducer of 55 ± 1 mm.

Flattener

Flattener + QHY600M with Filter Wheel and Off-Axis Guider



Placeholder image (must be replaced)


The Flattener is also from TS-Optics and maintains the focal length unchanged:

- Camera side M48x0.75 external thread with 55mm working distance
- Camera side M54x0.75 internal thread with 62.5mm working distance
- Camera side M78x1 internal thread with 77.5mm working distance
- Telescope side M92x1 external thread

Position	Adapter	Optical Path Length [mm]	
1	TS Optics M78a → M68a Adapter	3.3	
2	M68a → M68i Conversion Ring	12	
3	M68 Fine-Tuning Ring (golden)	1	
4	M68 Fine-Tuning Ring (black), optional	0.3	Adapter combination as in the table (ascending from left to right)
5	M68a → M62a Adapter	5	

As mentioned earlier, the working distance from the M78 thread is 77.5mm, with the camera along with the filter wheel, filter, and off-axis guider having a back focus of 56mm. The remaining 21.5mm are bridged by the above adapter combination. Last update: 2024/02/19 en:ost:telescope:ts-apo https://polaris.astro.physik.uni-potsdam.de/wiki/doku.php?id=en:ost:telescope:ts-apo&rev=1708382738 22:45

Flattener + QHY600M with UFC Filter Holder

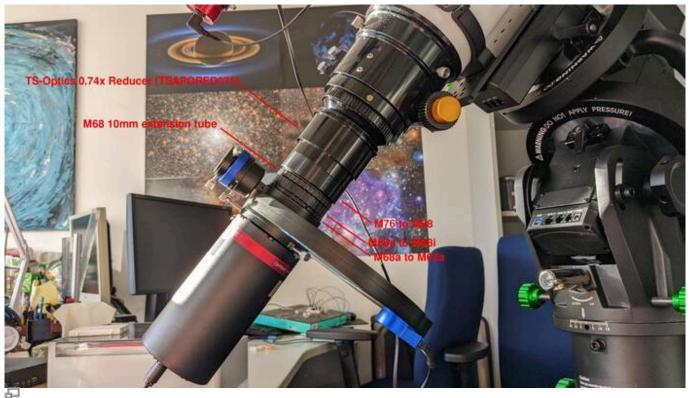

Placeholder image (must be replaced)

Position	Adapter	Optical Path Length [mm]	
1	TS Optics M78a → M68a Adapter	3.3	
2	M68 Fine-Tuning Ring (golden)	1	
3	M68 10mm Extension	10	
4	M68 Fine-Tuning Ring (golden)	1	Placeholder image (to be replaced)
5	M68 Fine-Tuning Ring (red), optional	0.5	
n I	M68a → M68i Conversion Ring	12	
7	M68 Fine-Tuning Ring (golden)	1	
8	M68 Fine-Tuning Ring (black), optional	0.3	
9	UFC S70 / M68x1	1	
10	UFC Tilter	9.75-10.5	
11	UFC Base	13]
12	Camera Adapter M54 → Camera Clamp	6	

Together, this results in an optical path length of approximately 58.8 mm depending on the tilter

position and installed fine-tuning rings. Combined with the 17.5 mm back focus of the QHY600M and the installed filter, this leads to 77.5 mm from the M78 thread.

Solar Observations

Solar Filter

A tension-free mounted solar filter for visual observations from Baader Planetarium is used. This filter is simply mounted in front of the objective. Any eyepiece and camera can then be used. However, the Herschel prism described below usually allows for better imaging quality.



APO with Herschel Prism

Herschel Prism

We also have a Herschel Prism Mark II from Baader Planetarium, which allows us to observe the sun in white light with excellent quality. The Herschel prism can be equipped with different filters, allowing it to be configured for visual observations as well as camera observations with very short exposure times. Details about the Herschel prism and its operation can be found in the associated article.

The Herschel prism can be used in many different configurations, for example, with and without a Barlow lens. However, the standard procedure is to insert it directly into the 2" eyepiece clamp of the APO.



Herschel Prism Mark II

APO with Herschel Prism

Last update: 2024/02/19 en:ost:telescope:ts-apo.https://polaris.astro.physik.uni-potsdam.de/wiki/doku.php?id=en:ost:telescope:ts-apo&rev=1708382738 22:45

Different models can also be used on the camera side. The prism has a 2" eyepiece clamp for holding corresponding accessories.

From: https://polaris.astro.physik.uni-potsdam.de/wiki/ - OST Wiki

Permanent link: https://polaris.astro.physik.uni-potsdam.de/wiki/doku.php?id=en:ost:telescope:ts-apo&rev=1708382738

Last update: 2024/02/19 22:45

